

Fitting a Replacement Filter collar

This entails the use of acetone in a syringe. Normal precautions should be taken including wearing protective glasses.

1. Remove the screw on filter from the collar – this is not always the case in these pictures.
2. Peel back the orange collar and insert the tie the correct way round. That is ensuring that when returned to the correct position the tie can be pulled tight. Do not tighten the tie.



Make sure the collar will fit over the precision mesh. It is better to rehearse fitting over the mesh, folding the orange rim back down over the collar and being ready to tighten the tie. When the acetone is applied there are only 5 seconds before the collar has to be in place.

3. Hold the collar away from you and apply acetone from the syringe all around the edge of the collar. Do not flood the collar or use excess amounts so that it drips.
4. Immediately insert the precision mesh inside the collar, reverse the orange material and pull down as far as possible.
5. Pull the tie very tight once you are sure there are no creases in the orange material and the precision mesh is in the right place.

6. Cut off the excess from the tie being very careful to avoid sharp edges on the net material.
7. Refit the screw on filter and tighten.